Clearing Biohazard Waste

Those on duty:

- 1. Clear Small bins on tables and bins in hood. If bags are wet and broken, please replace.
- 2. Clear waste from bin near sink into large bin near bacterial hood
- 3. Double bagging is required IF bags are broken/leaking
- 4. When bags are 70-80% full, tape the mouth shut and place in large biohazard disposal bin
- 5. Wash up and reward yourself! =)

Not on duty:

- 1. Clear Small bins if overflowing
- 2. Clear bins in hoods immediately after use
- 3. Decontaminate liquid biological waste using Presept tablets
- 4. We have no cytotoxic waste. If Cytotoxic waste is created, please contact Yi Han or Ivesc with explanation
- 5. If biological spills occur, refer to SOPs on biological waste management and use the biological spill kit under the sink. Please contact the relevant authorities (on main door)
- 6. Thank you for helping to keep the lab safe for all!

